

# Corgel™ BioHydrogel (TS-NaHy)

## Corgel™ Cross-linking Procedure

### CAUTION:

**PLEASE READ ALL INSTRUCTIONS PRIOR TO USE,  
AS THIS GEL SETS UP QUICKLY.**

### Product Description/Chemistry

Corgel™ tyramine substituted sodium hyaluronate (TS-NaHy), is a novel hyaluronan-based hydrogel system formed by either *in vitro* or *in situ* cross-linking driven by peroxidase.



**Illustration of the Cross-linking of TS-NaHy - not reflective of the cross-linking density**

*The use of this protocol will require optimization by the end user.*

**PLEASE REFRIGERATE UPON ARRIVAL**

### Reagents Provided

- Tyramine-Substituted Sodium Hyaluronate (TS-NaHy) (X% Substitution - **see bottle label**) - 250 mg powder. (Storage: 2 - 8° C)
- Horseradish Peroxidase (HRP) in PBS (10 U/mL) - 25 mL. Avoid prolonged exposure to light. (Storage: 2 - 8°C)
- Hydrogen Peroxide (H<sub>2</sub>O<sub>2</sub>) in PBS (1.0%) - 5 mL. Avoid prolonged exposure to light. (Storage: 2 - 8°C)

### Expiration

- Lifecore has not established an expiration date for the product.

### Procedures

*All steps may be performed at ambient temperature, unless otherwise specified.*

**Caution** - If you require more than one concentration of hydrogel, either aliquot the TS-NaHy before hydration or hydrate to the highest desired concentration and then aliquot to dilute to the desired concentrations.

1. Add HRP solution to hydrate TS-NaHy to the desired concentration. Use sufficient agitation and/or time to ensure that the solution is fully hydrated. When targeting a final gel concentration of TS-NaHy, take into account the amount of peroxide solution (step 2) to be added. Avoid direct sunlight.

**Caution** - Aliquot the TS-NaHy solution for dilution and/or make additions (for example, adding cells or drugs) as required before proceeding to the next step.

**Notes:**

- Gel concentrations below the lowest values and above the highest values in Table 1 may be generated, but the final hydrogel concentrations may not correlate due to expansion or contraction of the gel matrix.
  - The gel consistency will vary from a soft gel to firm gel particulates, and from optically clear to slightly colored gel, depending on TS-NaHy substitution, gel concentration, H<sub>2</sub>O<sub>2</sub> concentration, and additional components added.
  - The gel will set up within seconds, however maximum strength will be achieved after 16 hours.
  - Cells, bioactive agents, or other components may be added to the TS-NaHy solution during hydration, prior to H<sub>2</sub>O<sub>2</sub> addition, or after gel formation. These additives may affect the physical and biological properties of the gel.
2. Crosslink the TS-NaHy solution by adding H<sub>2</sub>O<sub>2</sub> solution of a desired concentration. The recommended volume ratio of H<sub>2</sub>O<sub>2</sub> solution to TS-NaHy solution is 1 to 25. Mix during addition if required. The H<sub>2</sub>O<sub>2</sub> reagent should be diluted with water to the desired concentration for cross-linking. Examples of the recommended H<sub>2</sub>O<sub>2</sub> concentrations are given in Table 1. Manual mixing may be applied to mix two components.

**Caution** - Use H<sub>2</sub>O<sub>2</sub> at a concentration not exceeding your predetermined limits.

**Table 1.** Examples of H<sub>2</sub>O<sub>2</sub> Concentration Required at Various Substitution percentages and Gel Concentrations

Substitution (%)	1.5	1.5	1.5
Gel concentration (mg/mL)	10	25	40
H <sub>2</sub> O <sub>2</sub> for cross-linking (%)	0.036	0.090	0.144

Substitution (%)	2.8	2.8	2.8
Gel concentration (mg/mL)	20	30	40
H <sub>2</sub> O <sub>2</sub> for cross-linking (%)	0.134	0.201	0.267

Substitution (%)	5.5	5.5	5.5
Gel concentration (mg/mL)	30	40	50
H <sub>2</sub> O <sub>2</sub> for cross-linking (%)	0.391	0.522	0.652

**Storage**

- Upon hydration, if not immediately used, TS-NaHy solution should be stored at: 2 – 8°C to minimize microbial growth. Prolonged storage is not recommended.
- If not immediately used, the crosslinked gel should be stored at 2 – 8°C to minimize microbial growth.

**FOR RESEARCH USE ONLY - NOT FOR HUMAN USE.**

Technology covered by U.S. PATENTS – 6,982,298; 7,368,502; 7,465,766 and related foreign patents.

**LIFECORE**

BIOMEDICAL Lifecore Biomedical, LLC 3515 Lyman Boulevard, Chaska MN USA 55318  
www.lifecore.com 1.800.348.4368 / 952.368.6321